









# 冷房安全裝置操作說明

## Instructions for Using Safety Devices in the Cold Room

1. 冷房門不會自動上鎖，必須從外面用力蓋上才會鎖上。  
The door of the cold room won't close automatically, so you must close it from outside.
2. 冷房門的內側設置安全推桿(彈簧門把)，由內往外推,門即開啟。  
Push the safety push-rod from inside to open the door.
3. 若安全推桿失效, 請立即逆時針(向左)旋轉緊急逃生鈕到底, 讓鎖頭鬆脫, 由內往外推, 門即開啟;或按下緊急警報鈕，警鈴作響燈號警示，等候救援。  
If the safety push-rod fails:
  - Turn the emergency door opening device to the left, and then push the door, or
  - Press the emergency button

1. 非自動上鎖	2.安全推桿	3. 緊急逃生鈕	4. 緊急警報鈕
 <p>非自動上鎖</p>	 <p>安全推桿由內向外推</p>	 <p>緊急逃生鈕左旋到底</p>	 <p>按下警報鈕</p>
The door won't lock automatically.	Push the safety push-rod	Turn the emergency door opening device to the left.	Press the emergency button.
		  <p>外側鎖頭鬆脫</p>	  <p>警鈴作響 燈號警示</p>
		To unlock the lock.	The alarm and signal will be turned on.